

***CETP* Activity Assay**

REF ELI-6279

1. Introduction

Cholesteryl ester transfer protein (CETP) is a member of the lipid transfer/lipoplysaccharide binding protein gene family. CETP transfers neutral lipids from high-density lipoprotein (HDL) to very low-density lipoprotein (VLDL) and is present in normal human plasma and serum.

The CETP Activity Assay Kit uses a donor molecule containing a fluorescent self-quenched neutral lipid that is transferred to an acceptor molecule in the presence of CETP. CETP-mediated transfer of the fluorescent neutral lipid to the acceptor molecule results in an increase in fluorescence (Excitation: 465 nm; Emission: 535 nm).

2. Reagents and Materials Supplied

100 assays; Store at 4°C

Component	Volume	Cap color
Donor Molecule	1 ml	Green
Acceptor Molecule	1 ml	Blue
CETP Assay Buffer (10X)	5 ml	Clear
Positive Control (Rabbit Serum)	30 µl	Red

3. Procedure of the Test

I. General Consideration for Using Fluorometer and Plate Reader

We recommend using a microtiter plate for the assay. The microtiter plates should be sealed as tightly as possible with plate sealer and incubated in a sealed, humidified chamber to prevent evaporation.

If using a regular fluorometer for sample reading, the samples should be diluted to 500 µl with 1X CETP Assay Buffer before read.

II. Preparation of Standard Curve

- Standard curve is prepared by making serial dilutions of the donor molecule in isopropanol and subsequently recording the fluorescence intensity of each dilution, using isopropanol alone as a blank.
- Prepare 6 test tubes labeled T0 to T5, each contains 0.2 ml of isopropanol; the tube labeled T5 should contain an additional 0.2 ml of isopropanol.
- Add 2 µl Donor Molecule to T5, vortex to mix well.
- Transfer 0.2 ml from T5 to T4. Mix and then transfer 0.2 ml from T4 to T3. Mix and then transfer 0.2 ml from T3 to T2. Mix and then transfer 0.2 ml from T2 to T1. The Donor Molecule solution contains 0.1 mM labeled lipids and thus the standard curve samples contain 0, 6.25, 12.5, 25, 50, 100 pmol donor molecule.
- Read the fluorescence intensity (Ex. = 465 nm; Em. = 535 nm) of the samples from T0 to T5.
- Apply the fluorescence intensity values of the standard curve directly to your results to express activity of the plasma sample (pmoles/ml plasma/hour), or use Relative Fluorescence Units (RFU)/µl plasma/hour.

Assay Procedure

- For each reaction, add the following components:
 - 10 µl Donor Molecule
 - 10 µl Acceptor Molecule
 - 20 µl 10X CETP Assay Buffer

1-3 µl Your Sample (serum or plasma)
ddH₂O To a total of 200 µl

For positive control, add 1-3 µl of Rabbit Serum instead of your sample.

Prepare a blank that contains no CETP Source.

- Incubate for 30-60 minutes at 37°C.
- Measure the fluorescence intensity of the blank, samples, and positive control using a fluorescence plate reader or fluorometer (Ex. = 465 nm; Em. = 535 nm). Due to the nature of the self-quenched probe, background fluorescence can be significant; therefore, fluorescence intensity from each sample should be corrected by subtracting the blank fluorescence intensity. The increase in fluorescence intensity is usually 0.2-2 fold over blank.
- Calculate the activity of the plasma sample:

$$Y = MX + B$$


Where:

Y = Fluorescence Intensity of Sample
– Fluorescence Intensity of Blank
M = Slope of the Standard Curve
X = Concentration of Plasma Sample
B = Intercept

Example:

Y = 16283 – 8655 = 7628
M = 74.5
B = 480
7628 = 74.5X + 480
X = 95.9 pmol/µl sample/1 hour

CETP Activity Assay with

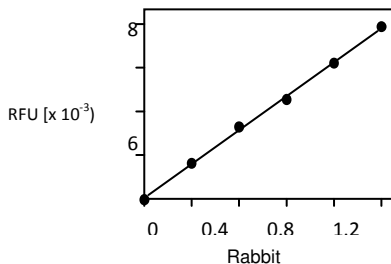











4. Troubleshooting

Problems	Cause	Solution
Assay not working	Use of ice-cold assay buffer	Assay buffer must be at room temperature
	Omission of a step in the protocol	Refer and follow the data sheet precisely
	Plate read at incorrect wavelength	Check the wavelength in the data sheet and the filter settings of the instrument
	Use of a different 96-well plate	Fluorescence: Black plates (clear bottoms) Luminescence: White plates; Colorimeters: Clear plates
Samples with erratic readings	Use of an incompatible sample type	Refer to data sheet for details about incompatible samples
	Samples prepared in a different buffer	Use the assay buffer provided in the kit or refer to data sheet for instructions
	Cell/ tissue samples were not completely homogenized	Use Dounce homogenizer (increase the number of strokes) observe for lysis under microscope
	Samples used after multiple free-thaw cycles	Aliquot and freeze samples if needed to use multiple times
	Presence of interfering substance in the sample	Troubleshoot if needed
	Use of old or inappropriately stored samples	Use fresh samples or store at correct temperatures until use
Lower/ Higher readings in Samples and Standards	Improperly thawed components	Thaw all components completely and mix gently before use
	Use of expired kit or improperly stored reagents	Always check the expiry date and store the components appropriately
	Allowing the reagents to sit for extended times on ice	Always thaw and prepare fresh reaction mix before use
	Incorrect incubation times or temperatures	Refer to datasheet & verify correct incubation times and temperatures
	Incorrect volumes used	Use calibrated pipettes and aliquot correctly
Readings do not follow a linear pattern for Standard curve	Use of partially thawed components	Thaw and resuspend all components before preparing the reaction mix
	Pipetting errors in the standard	Avoid pipetting small volumes
	Pipetting errors	Prepare a master reaction mix

	in the reaction mix	whenever possible
	Air bubbles formed in well	Pipette gently against the wall of the tubes
	Standard stock is at an incorrect concentration	Always refer to the dilutions in the data sheet
	Calculation errors	Recheck calculations after referring to the data sheet
	Substituting reagents from older kits / lots	Use fresh components from the same kit
Unanticipated results	Measured at incorrect wavelength	Check the equipment and the filter setting
	Samples contain interfering substances	Troubleshoot if it interferes with the kit
	Use of incompatible sample type	Refer to data sheet to check if sample is compatible with the kit or optimization is needed
	Sample readings above/below the linear range	Concentrate / Dilute sample so as to be in the linear range

Rev. 1, 2012-10-24 SS

Symbol	Deutsch	English	Français	Español	Italiano	Polski
	CE Konformitätszeichen	CE marking of conformity	Conformité aux normes européennes	Conformidad europea	Conformità europea	Znak zgodności CE
	Gebrauchsanweisung beachten	Consult instructions for use	Consulter la notice d'utilisation	Consultense las instrucciones de uso	Consultare le istruzioni per l'uso	Przestrzegać instrukcji obsługi
	In-vitro-Diagnostika	In-vitro diagnostic medical device	Dispositif médical de diagnostic in vitro	Producto sanitario para diagnóstico in vitro	Dispositivo medico-diagnostico in vitro	Tylko do diagnostyki in vitro
	Temperaturbegrenzung	Temperature limitation	Limites de température	Limitación de temperatura	Limiti di temperatura	Temperatura przechowywania
	Chargenbezeichnung	Batch code	Code du lot	Código de lote	Codice lotto	Numer serii
	Nicht zur Wiederverwendung	Do not reuse	Ne pas réutiliser	No reutilizar	Non riutilizzare	Tylko do jednorazowego użytku
	Verwendbar bis	Use by	Utiliser jusqu'au	Fecha de caducidad	Utilizzare entro	Data ważności
	Bestellnummer	Catalogue Number	Référence du catalogue	Número de catálogo	Riferimento di Catalogo	Numer katalogowy
	Hersteller	Manufacturer	Fabricant	Fabricante	Fabbricante	Producent
	Ausreichend für <n> Ansätze	Sufficient for <n> tests	Suffisant pour pour "n" tests	Válido para para <n> ensayos	Sufficiente per "n" saggi	Wystarczający na <n> Powtórzeń

Our Teams

Germany:

Regensburg

Tel: +49 941 290 10-0
 Fax: +49 941 290 10-50

Moers

Tel: +49 2841 99820-0
 Fax: +49 2841 99820-1

Austria:

Tel: +49 941 290 10-29
 Free Tel: 0800 291 565
 Fax: +49 290 10-50
 Free Fax: 0800 298 197

UK & Ireland:

Tel: +49 941 290 10-18
 Free Tel – UK: 0800 234 1237
 Free Tel – IRE: 1800 555 080
 Fax: +49 290 10-50

France:

France Tel: 0800 915 240
 France Fax: 0800 909 493

Switzerland

Swiss Tel: 0800 564 720
 Swiss Fax: 0800 837 476

Belgium

Belgium Tel: 0800 718 82
 Belgium Fax: 0800 747 07

Luxembourg

Lux. Tel: 800 211 16
 Lux. Fax: 800 261 79

Spain:

Tel: +49 941 290 10-759
 Free Tel: 900 938 315
 Fax: +49 941 290 10-50
 Free Fax: 900 984 992

Italy:

Tel: +49 941 290 10-34
 Fax: +49 941 290 10-50

Poland:

Tel: +49 941 290 10-44
 Free Tel: 00 800 491 15 95
 Fax: +49 941 290 10-50
 Free Fax: 00 800 491 15 94

Netherlands:

Tel: +31 30 75 600
 Free Tel: 0800 0222 890
 Fax: +31 70 30 30 775
 Free Fax: 0800 024 9519

Denmark:

Tel: +31 703075 603
 Free Tel: +45 80 88 87 53
 Tax: +31 703030 775

Laboratory Diagnostics Team:

Tel: +49 941 290 10-40
 Fax: +49 941 290 10-50



nal von minden GmbH
 Carl-Zeiss-Strasse 12
 47445 Moers
 Germany

www.nal-vonminden.com
 info@nal-vonminden.com
 Fon: +49 2841 99820-0
 Fax: +49 2841 99820-1