

Caspase-Family Fluorometric Substrate Set

CATALOG #: K131-7-25

LOT #: _____

STORAGE CONDITIONS: Store at -20°C.

SHELF LIFE: 1 year under proper storage conditions

DESCRIPTION:

Fluorometric substrates for assaying activities of members of caspase family proteases. All substrates are provided in liquid ready-to-use form.

Kit CONTENTS: Each of the following substrates is dissolved in DMSO

Concentration	Description	Volume	Part Number
1 mM	Caspase-1 Substrate, Ac-YVAD-AFC	125 µl	K110-25-3
1 mM	Caspase-2 Substrate, Ac-VDVAD-AFC	125 µl	K116-25-3
1 mM	Caspase-3 Substrate, Ac-DEVD-AFC	125 µl	K105-25-3
1 mM	Caspase-5 Substrate, Ac-WEHD-AFC	125 µl	K122-25-3
1 mM	Caspase-6 Substrate, Ac-VEID-AFC	125 µl	K114-25-3
1 mM	Caspase-8 Substrate, Ac-IETD-AFC	125 µl	K112-25-3
1 mM	Caspase-9 Substrate, Ac-LEHD-AFC	125 µl	K118-25-3

ASSAY PROCEDURE:

1. Induce apoptosis in cells by desired method. Concurrently incubate a control culture *without* induction.
2. Count cells and pellet 1-5 x 10⁶ cells or use 50-200 µg cell lysates if protein concentration has been measured.
3. Resuspend cells in 50 µl of chilled Cell Lysis Buffer (Cat.# 1067-100).
4. Incubate cells on ice for 10 minutes.
5. Add 50 µl of 2X Reaction Buffer (Cat. # 1068-20) containing 10 mM DTT (Cat.# 1201-1) to each sample.
6. Add 5 µl of the 1 mM AFC conjugated substrates (50 µM final conc.) into each tube individually and incubate at 37°C for 1-2 hour.
7. Read samples in a fluorometer equipped with a 400-nm excitation filter and 505-nm emission filter. For a plate-reading set-up, transfer the samples to a 96-well plate. You may perform the entire assay directly in a 96-well plate.
Fold-increase in caspase activity can be determined by comparing these results with the level of the uninduced control.

FOR RESEARCH USE ONLY! Not to be used on humans.

RELATED PRODUCTS:

Apoptosis Detection Kits & Reagents

- Annexin V Kits & Bulk Reagents
- Caspase Assay Kits & Reagents
- Mitochondrial Apoptosis Kits & Reagents
- Nuclear Apoptosis Kits & Reagents
- Apoptosis Inducers and Set
- Apoptosis siRNA Vectors

Cell Fractionation System

- Mitochondria/Cytosol Fractionation Kit
- Nuclear/Cytosol Fractionation Kit
- Membrane Protein Extraction Kit
- Cytosol/Particulate Rapid Separation Kit
- Mammalian Cell Extraction Kit
- FractionPREP Fractionation System

Cell Proliferation & Senescence

- Quick Cell Proliferation Assay Kit
- Senescence Detection Kit
- High Throughput Apoptosis/Cell Viability Assay Kits
- LDH-Cytotoxicity Assay Kit
- Bioluminescence Cytotoxicity Assay Kit
- Live/Dead Cell Staining Kit

Cell Damage & Repair

- HDAC Fluorometric & Colorimetric Assays & Drug Discovery Kits
- HAT Colorimetric Assay Kit & Reagents
- DNA Damage Quantification Kit
- Glutathione & Nitric Oxide Fluorometric & Colorimetric Assay Kits

Signal Transduction

- cAMP & cGMP Assay Kits
- Akt & JNK Activity Assay Kits
- Beta-Secretase Activity Assay Kit

Adipocyte & Lipid Transfer

- Recombinant Adiponectin, Survivin, & Leptin
- CETP Activity Assay & Drug Discovery Kits
- PLTP Activity Assay & Drug Discovery Kits
- Total Cholesterol Quantification Kit

Molecular Biology & Reporter Assays

- siRNA Vectors
- Cloning Insert Quick Screening Kit
- Mitochondrial & Genomic DNA Isolation Kits
- 5 Minutes DNA Ligation Kit
- 20 Minutes Gel Staining/Destaining Kit
- β -Galactosidase Staining Kit & Luciferase Reporter Assay Kit

Growth Factors and Cytokines

GENERAL TROUBLESHOOTING GUIDE FOR CASPASE COLORIMETRIC AND FLUOROMETRIC KITS:

Problems	Cause	Solution
Assay not working	<ul style="list-style-type: none"> • Cells did not lyse completely • Experiment was not performed at optimal time after apoptosis induction • Plate read at incorrect wavelength • Old DTT used 	<ul style="list-style-type: none"> • Resuspend the cell pellet in the lysis buffer and incubate as described in the datasheet • Perform a time-course induction experiment for apoptosis • Check the wavelength listed in the datasheet and the filter settings of the instrument • Always use freshly thawed DTT in the cell lysis buffer
High Background	<ul style="list-style-type: none"> • Increased amount of cell lysate used • Increased amounts of components added due to incorrect pipetting • Incubation of cell samples for extended periods • Use of expired kit or improperly stored reagents • Contaminated cells 	<ul style="list-style-type: none"> • Refer to datasheet and use the suggested cell number to prepare lysates • Use calibrated pipettes • Refer to datasheet and incubate for exact times • Always check the expiry date and store the individual components appropriately • Check for bacterial/ yeast/ mycoplasma contamination
Lower signal levels	<ul style="list-style-type: none"> • Cells did not initiate apoptosis • Very few cells used for analysis • Use of samples stored for a long time • Incorrect setting of the equipment used to read samples • Allowing the reagents to sit for extended times on ice 	<ul style="list-style-type: none"> • Determine the time-point for initiation of apoptosis after induction (time-course experiment) • Refer to datasheet for appropriate cell number • Use fresh samples or aliquot and store and use within one month for the assay • Refer to datasheet and use the recommended filter setting • Always thaw and prepare fresh reaction mix before use
Samples with erratic readings	<ul style="list-style-type: none"> • Uneven number of cells seeded in the wells • Samples prepared in a different buffer • Adherent cells dislodged and lost at the time of experiment • Cell/ tissue samples were not completely homogenized • Samples used after multiple freeze-thaw cycles • Presence of interfering substance in the sample • Use of old or inappropriately stored samples 	<ul style="list-style-type: none"> • Seed only equal number of healthy cells (correct passage number) • Use the cell lysis buffer provided in the kit • Perform experiment gently and in duplicates/triplicates; apoptotic cells may become floaters • Use Dounce homogenizer (increase the number of strokes); observe efficiency of lysis under microscope • Aliquot and freeze samples, if needed to use multiple times • Troubleshoot as needed • Use fresh samples or store at correct temperatures until use
Unanticipated results	<ul style="list-style-type: none"> • Measured at incorrect wavelength • Cell samples contain interfering substances 	<ul style="list-style-type: none"> • Check the equipment and the filter setting • Troubleshoot if it interferes with the kit (run proper controls)
General issues	<ul style="list-style-type: none"> • Improperly thawed components • Incorrect incubation times or temperatures • Incorrect volumes used • Air bubbles formed in the well/tube • Substituting reagents from older kits/ lots • Use of a different 96-well plate 	<ul style="list-style-type: none"> • Thaw all components completely and mix gently before use • Refer to datasheet & verify the correct incubation times and temperatures • Use calibrated pipettes and aliquot correctly • Pipette gently against the wall of the well/tubes • Use fresh components from the same kit • Fluorescence: Black plates; Absorbance: Clear plates
Note: The most probable cause is listed under each section. Causes may overlap with other sections.		