


New 8-OHdG Check ELISA

ELISA kit for 8-hydroxy-2'-deoxyguanosine

FOR RESEARCH USE ONLY

About 8-hydroxy-2'-deoxyguanosine (8-OHdG):

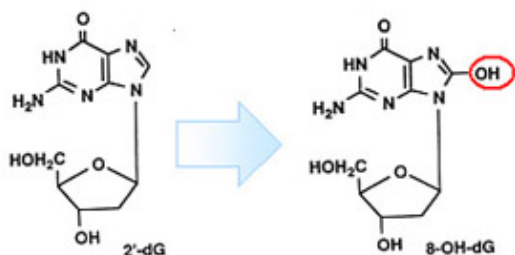

Formation of 8-hydroxy-2'-deoxyguanosine (8-OHdG) by oxygen radicals.
H Kasai: Environmental Mutagen Research, Vol. 10, p73-78 (1988)

8-hydroxy-2'-deoxyguanosine (8-OHdG) is a product of oxidatively damaged DNA formed by hydroxy radical, singlet oxygen and direct photodynamic action. 8-OHdG can be detected in tissue, serum, urine and other biomaterials. New 8-OHdG Check is a competitive enzyme-linked immunosorbent assay (ELISA) utilising monoclonal antibody (clone N45.1) which is highly specific for DNA damage, not cross react with RNA oxidation products such as 8-hydroxy-guanine and 8-hydroxy-guanosine. This product is suitable for detection of 8-OHdG in urine and other biomaterials from human and animals.



This product is a 8-OHdG ELISA kit utilizing anti 8-OHdG monoclonal antibody (clone N45.1) which is highly specific for 8-OHdG. We provide two types of 8-OHdG ELISA kits with different assay range. New 8-OHdG Check ELISA is suitable for urine and serum sample from animal and human. If you are planning to measure 8-OHdG in human serum, tissue, cultured cells, we recommend to use 'Highly Sensitive 8-OHdG Check ELISA'.

Specifications

Assay principle: Competitive ELISA (detection: 450 nm)

Specificity: Specific for 8-OHdG. Antibody have been tested to 8-OHdG analogues (guanosine(G), 7-methyl-G, 6-SH-G, 8-Bromo-G, dA, dC, dT, dI, dU, dG, O6-methyl-dG, 8-OHdA, guanine(Gua), O6-methyl-Gua, 8-OH-Gua, uric acid, urea, creatine, creatinine, 8-sulfhydryl-G, 8-OH-G).

Measuring range: 0.5 to 200 ng/mL

Format: 96 wells (18 samples in triple assays)

Applications: Urine from human and animals.
For serum and tissue samples, Highly Sensitive 8-OHdG Check (code #KOG-HS10E) is recommended.

Storage: Store at 4 - 10°C (don't freeze).

Product name	Code	Assay range	Format	Application
New 8-OHdG Check	KOG-200SE	0.5-200 ng/mL	96 wells	Urine and animal serum.
New 8-OHdG Check Trial Package	KOG-200TE	0.5-200 ng/mL	32 wells	

Related products: [Highly sensitive 8-OHdG Check ELISA](#) [Anti 8-OHdG monoclonal antibody](#)

FOR RESEARCH USE ONLY. Not for diagnostic, medical or other use.

Japan Institute for the Control of Aging (JaiCA) TEL:+81-538490125 / FAX:+81-538491267 <http://www.jaica.com>

Gentaur Molecular Products
Voortstraat 49
1910 Kampenhout, Belgium

Content of this kit

8-OHdG Microtiter Plate:	Precoated with 8-OHdG(12 X 8w ells, split type)*	1 plate
Primary Antibody:	Anti 8-OHdG antibody, pow der.	1 vial
Primary Antibody Solution		1 vial (6mL)
Secondary Antibody:	HRP-anti mouse antibody, pow der.	1 vial
Secondary Antibody Solution:		1 vial (12mL)
Chromatic Solution:	3,3',5,5'-tetramethylbenzidine	1 vial (0.25mL)
Diluting Solution:	H ₂ O ₂ containing buffer.	1 vial (12mL)
Washing Solution(5x):		2 vials (26mL x 2)*
Reaction Terminating Solution:	1M Phosphoric acid.	1 vial (12mL)
Standard 8-OHdG Solution:	Purified 8-OHdG (0.5, 2, 8, 20, 80, 200 ng/mL)	1 vial each
Plate Seal:		2 sheets

*:Trial package (KOG-200TE) contains 32 w ells (4 X 8 w ells), and one vial of Washing solution(5X).

References

- 1) H.Kasai, P.F.Crain, Y.Kuchino, S.Nishimura,A.Ootsuyama and H.Tanooka : Formation of 8-hydroxyguanine moiety in cellular DNA by agents producing oxygen radicals and evidence for its repair. *Carcinogenesis* 7(11), p1849-1851 (1986)
[[8-OHdG/8-OHG is formed by reactive oxygen radicals. \]](#)
- 2) R.G.Cutler: Antioxidants and aging. *Am J Clin Nutr* 50,p373S-379S (1991)
[[Relationship between lifespan and antioxidants, enzymes and 8-OHdG. \]](#)
- 3) S.Toyokuni, T.Tanaka, Y.Hattori, Y.Nishiyama, A.Yoshida, K.Uchida, H.Hiai, H.Ochi and T.Osawa : Quantitative immunohistochemical determination of 8-hydroxy-2'-deoxyguanosine by a monoclonal antibody N45.1: Its application to ferric nitrilotriacetate-induced renal carcinogenesis model. *Lab Invest* 76(3), p365-374 (1997)
[[Specifity of the antibody clone N45.1. \]](#)
- 4) Saito S, Yamauchi H, Hasui Y, Kurashige J, Ochi H, Yoshida K: Quantitative determination of urinary 8-hydroxydeoxyguanosine (8-OH-dg) by using ELISA. *Res Commun Mol Pathol Pharmacol* 107(1-2),p39-44 (2000)
[[Development and assessment of 8-OHdG Check ELISA. \]](#)
- 5) Nagayoshi Y, Kawano H, Hokamaki J, Miyamoto S, Kojima S, Shimomura H, Tsujita K, Sakamoto T, Yoshimura M, Ogawa H: Urinary 8-hydroxy-2'-deoxyguanosine levels increase after reperfusion in acute myocardial infarction and may predict subsequent cardiac events. *Am J Cardiol*. 95(4), p514-517 (2005)
[[Measurement of 8-OHdG in human urine. \]](#)
- 6) Morillas-Ruiz J, Zafrilla P, Almar M, Cuevas MJ, Lopez FJ, Abellan P, Villegas JA, Gonzalez-Gallego J: The effects of an antioxidant-supplemented beverage on exercise-induced oxidative stress: results from a placebo-controlled double-blind study in cyclists. *Eur J Appl Physiol* 95(5-6),p543-549(2005)
[[Study about exercise, antioxidant supplements and urinary 8-OHdG. \]](#)
- 7) Someya T, Kaneko K, Yamada T, Yamashiro Y: Effect of a novel free radical scavenger, edaravone, on puromycin aminonucleoside induced nephrosis in rats. *Pediatr Nephrol* 20(10),p1430-1434(2005)
[[Effect of antioxidants to 8-OHdG in rat urine. \]](#)
- 8) Shiihara T, Kato M, Ichiyama T, Takahashi Y, Tanuma N, Miyata R, Hayasaka K: Acute encephalopathy with refractory status epilepticus: Bilateral mesial temporal and claustral lesions, associated with a peripheral marker of oxidative DNA damage. *J Neurol Sci* 250(1-2),p159-161(2006)
[[Measurement of 8-OHdG in human urine, serum and cerebro spinal fluid\(CSF\). \]](#)

FOR RESEARCH USE ONLY. Not for diagnostic, medical or other use.

